



## Product Information

Trypsin (0.25 %) in DPBS (1x)

Catalog no: GBT01/01F

### General Information

Trypsin solutions are utilized to separate cells that are adhered to culture surfaces. They contain naturally occurring swine pancreatic trypsin. The amount of trypsin necessary to separate cells from their substrate is most strongly influenced by the kind of cell and the age of the culture. Numerous formulations need to be assessed in order to identify the best product for a certain application.

### Specification

Appearance	Clear frozen liquid
Storage and shelf life	Store at $\leq -15^{\circ}\text{C}$ . Avoid repeated freeze-thaw cycles. Preparation of aliquots recommended. Once opened, store at $4^{\circ}\text{C}$ and use within 2-4 weeks.
Shipping conditions	Frozen (Dry ice)
Thawing	$+37^{\circ}\text{C}$ water bath or overnight at $+2^{\circ}\text{C}$ to $+8^{\circ}\text{C}$ . Swirl gently to homogenize.

### Formulation

Components	Concentration mg/L
KCl	200.00
$\text{KH}_2\text{PO}_4$	200.00
NaCl	8000.00
$\text{Na}_2\text{HPO}_4$	1150.00
Trypsin	2500.00

### Guidelines for Use

#### Detachment of adherent cells:

Trypsin (0.25%) in DPBS (1x) solution is offered as a frozen, sterile liquid that is ready for use. The entire process should be carried out in a laminar flow hood with strict asepsis.

1. The product can be defrosted overnight at  $+2^{\circ}\text{C}$  to  $+8^{\circ}\text{C}$  or in a water bath at  $+37^{\circ}\text{C}$ .
2. Carefully aspirate the cell culture flask's whole media supply.
3. Rinse the cells in a salt solution devoid of  $\text{Ca}^{2+}$  and  $\text{Mg}^{2+}$  (see related products), aspirate, and discard.
4. Warm the trypsin solution up in a water bath at  $+37^{\circ}\text{C}$ . Make sure the cells are thoroughly covered in trypsin solution.
5. Keep the flask at  $+37^{\circ}\text{C}$  or, for more delicate cultures, at  $+2^{\circ}\text{C}$  to  $+8^{\circ}\text{C}$  or room temperature.

#### Disclaimer:

User must ensure suitability of the product(s) in their application prior to use. Products conform solely to the information contained in this and other related Products are not intended for human or animal diagnostic or therapeutic use but for laboratory, research or further manufacturing use only, unless otherwise specified. Statements contained herein should not be considered as a warranty of any kind, expressed or implied, and no liability is accepted for infringement of any patents.



## Product Information

Trypsin (0.25 %) in DPBS (1x)

Catalog no: GBT01/01F

6. Cells will appear spherical under a microscope after the trypsinization procedure is finished, and the flask's solution will look hazy. To prevent overexposure, check the flask often. Trypsin exposure should be limited in duration since it might harm cells.

The kind of cell, the culture's age, population density, serum content in the growth media, and the interval since the last subculture all affect how long it takes for cells to separate from the culture surface.


7. Use a trypsin inhibitor or a serum-containing medium to neutralise trypsin. The trypsin-containing supernatant should be discarded after gently centrifuging the cell solution.

8. Use new media to resuspend the cell pellet, then count or cultivate the cells as needed.

### Need Help?

---

If you have any further queries, please feel free to email our cell culture specialists at [info@genexisbiotech.com](mailto:info@genexisbiotech.com)

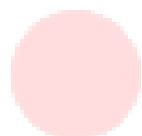

#### Disclaimer:

User must ensure suitability of the product(s) in their application prior to use. Products conform solely to the information contained in this and other related Products are not intended for human or animal diagnostic or therapeutic use but for laboratory, research or further manufacturing use only, unless otherwise specified. Statements contained herein should not be considered as a warranty of any kind, expressed or implied, and no liability is accepted for infringement of any patents.